



2F, No. 173, Shengli 6th St, Zhubei City, Hsinchu County 302, Taiwan  
Tel: +886 35721683 Fax: +886 35721633

# HRP Conjugation Kit

For simple and quick conjugation of IgG antibodies with HRP. For research use only - not intended for diagnostic use.

## Introduction:

HRP Conjugation Kit - provides a simple and rapid procedure to covalent labeling of antibodies with HRP by targeting primary amine groups (e.g. lysines). The kit simplifies the labeling process into a quick procedure, minimizing hands-on time. A key benefit is that the resulting conjugate is ready for use without needing post-conjugation purification, which helps maximize sample recovery and save time.

The antibody to be labeled should be purified, in an appropriate buffer for conjugation and at a suitable concentration conjugation (as described in buffer considerations).

The kit comes in 5 pack sizes:

- 3 x 10 µg containing 3 reactions (each for 10 - 20 µg)
- 100 µg containing 1 reaction (for 100 - 200 µg)
- 3 x 100 µg containing 3 reactions (each for 100 - 200 µg)
- 1 mg containing 1 reaction (for 1 - 2 mg)
- 3 x 1 mg containing 3 reactions (each for 1 - 2 mg)

Custom size conjugation kits up to 100 mg are available on demand. Please contact us to discuss your requirements.

## Materials Supplied:

Item	3 x 10 µg	100 µg	3 x 100 µg	1 mg	3 x 1 mg	Storage temperature
HRP mix	3 vials	1 vial	3 vials	1 vial	3 vials	-20°C
Modifier reagent	1 vial	1 vial	1 vial	1 vial	1 vial	-20°C
Quencher reagent	1 vial	1 vial	1 vial	1 vial	1 vial	-20°C

**Storage and Stability:** Store kit at -20 °C immediately upon receipt.

## Buffer considerations:

### Recommended pre-conjugation buffer components and conditions:

For optimal labeling efficiency and conjugate stability with the FITC Conjugation kit, the starting antibody or protein solution should ideally meet specific buffer requirements.

Recommended Buffer:

- Amine free buffer (e.g. MES, MOPS, HEPES, PBS, Borate buffer)
- pH:6.5-8.5
- Antibody concentration: 1-8mg/mL

**Note:** It is important to AVOID Nucleophilic components (primary amines e.g. amino acids or ethanolamine and thiols e.g. mercaptoethanol or DTT)

## Recommended starting amount and volume of antibody:

Vial Size	Lowest Amount of Antibody	Recommended Amount of Antibody (1:2 Antibody:HRP)	Maximum Amount of Antibody (1:1 Antibody:HRP)	Fixed Volume of Antibody
10 µg	Down to 10 µg at 1 mg/mL	Optimally 10 µg at 4 mg/mL	Up to 20 µg at 4 mg/mL	Fixed 10 µL
100 µg	Down to 100 µg at 1 mg/mL	Optimally 100 µg at 4 mg/mL	Up to 200 µg at 4 mg/mL	Fixed 100 µL
1 mg	Down to 1mg at 1 mg/mL	Optimally 1,000 µg at 4 mg/mL	Up to 2,000 µg at 4 mg/mL	Fixed 1,000 µL

### Note:

1. It is important to use the indicated fixed volumes. Modifying the reaction volume may result in poor quality conjugates.
2. Antibodies can be diluted using either milli-Q water or PBS.
3. Optimal conjugates are normally generated using the recommended amount of antibody, although using different amounts of antibody within the range provided will still generate quality conjugates.
4. Adding less than the recommended amount of antibody may result in unbound label post conjugation. This excess label will be deactivated by the quencher and removed during the first wash step of any application.

## Assay Procedure

Equilibrate all materials and prepared reagents to room temperature (20-25°C) prior to use.

### 1. Prepare the antibody:

Add 1 µL of Modifier reagent to each 10 µL of antibody to be labeled and mix gently.

### 2. Add antibody to the HRP:

Remove the cap from the vial containing the HRP Conjugation Mix. Pipette the prepared antibody sample (containing the Modifier reagent) directly onto the lyophilized material inside the vial. Resuspend gently by carefully withdrawing and re-dispensing the liquid once or twice using a pipette tip.

### 3. Incubate the reaction:

Replace the cap on the vial. Incubate the reaction mixture for either **3 hours at room temperature** (20-25°C) in the dark, or **overnight at 4°C**.

### 4. Quench the reaction:

After incubation, add 1 µL of Quencher reagent for every 10 µL of the initial antibody used, and mix gently. The conjugate will be ready for use after 30 minutes. The resulting conjugates do not require a purification step and can be used directly in downstream applications.

## Conjugate Storage

Typically, the antibody conjugate can be stored at 4°C for up to 12 months. For longer storage the conjugate can be stored at -20°C with a cryoprotectant such as 50% glycerol. The best storage conditions for any particular conjugate must be determined by experimentation and depend on the antibody stability and best storage conditions. The antibody conjugate should always be stored in the dark.