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Mouse Antibody Isotyping Kit Product Insert & Protocol

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- Purified Monoclonal Antibody: Dilute to 1 – 5 µg/mL.
- *(Note: Excessive concentration may cause a 'Hook Effect,' leading to faint or absent lines.)*

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IV. Assay Procedure

1. Equilibrate: Bring sealed pouch and buffer to room temperature (20°C - 30°C) for 15–30 minutes before testing.
2. Unseal: Open aluminum foil pouch and remove Test Cassettes A and B. Place them horizontally on a flat surface.
3. Load Sample: Pipette 80 – 100 µL of diluted sample vertically and slowly into the Sample Well of both Cassettes A and B.
4. Incubate: Leave undisturbed at room temperature for capillary migration.
5. Read Results: Read and record results within 5 – 10 minutes. Results after 15 minutes are invalid.

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V. Result Interpretation

- Valid Test: The Control Line (C) at the top of both Cassette A and Cassette B must show a distinct red band.
- Positive Result: If the C-line is normal, observe which specific isotype letter exhibits a red band. E.g., if G1 line on Cassette B appears, the isotype is IgG1.
- Multiple Bands (Mixed): If more than one isotype band appears simultaneously (e.g., both G1 and M), it indicates the hybridoma cell line may not be monoclonal yet or contaminated. Subcloning is recommended.
- Invalid Test: If no red band appears at the Control Line (C) on either cassette, the test has failed. Re-test with a new kit.

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VI. Precautions

1. For Research Use Only (RUO). Not for use in diagnostic procedures.
2. Test cassettes are moisture-sensitive. Use within 10 minutes after unsealing.
3. Do not mix or substitute components from different lot numbers.